## USB Driver installation procedure

Install two drivers.

- CP210x USB Composite Device
  - CP210x USB to UART Bridge Controller
- (Two drivers must be installed.)

"Win2K\_XP" folder will be copied in the hard disc from the CD-ROM.

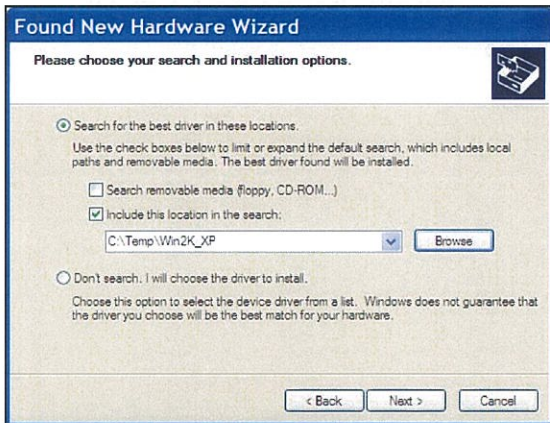
(In this explanation, copy to "C:\Temp")

Connect USB I/F(interface) cable to P.C.

The window of "Found New Hardware Wizard" will show up a short while.

Select "Install from a list or specific location". Click "Next".

Select "Search for the best driver...". Uncheck "Search removable media". Check "Include this location in the search". Click "Browse" and select "Win2K\_XP" folder. Click "OK". Click "Next".

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

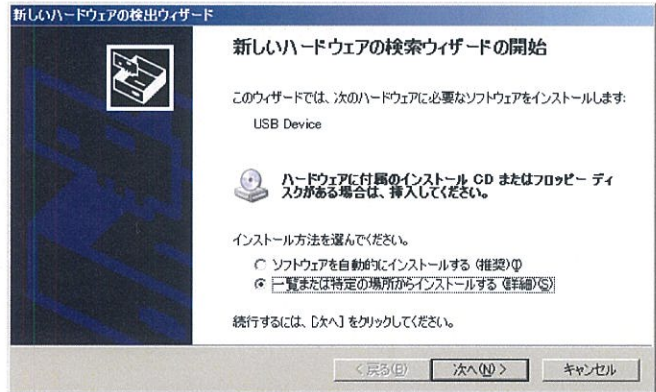
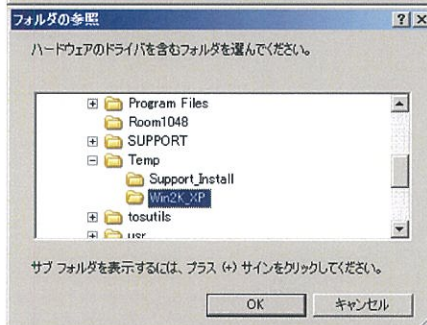
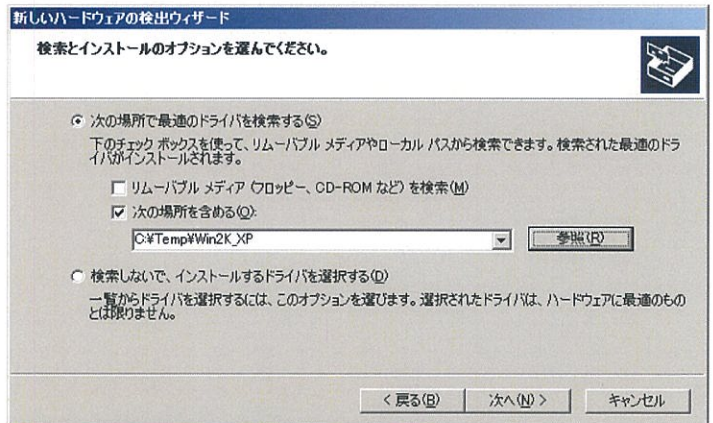
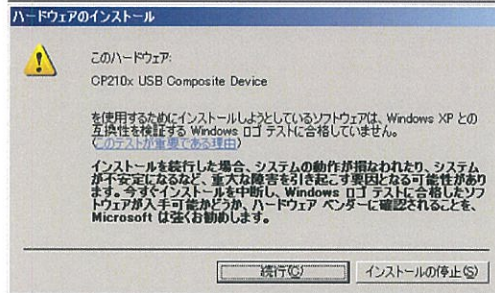
"CP210x USB Composite Device" installation begins. During the installation, the window would show up, click "Continue anyway".

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"CP210x USB Composite Device" installation finished. Click "Finish".

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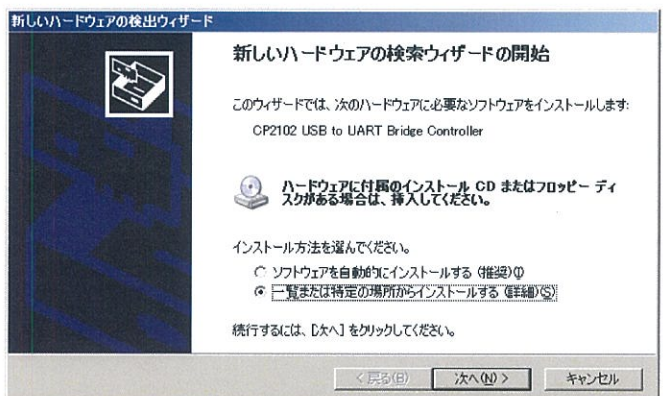
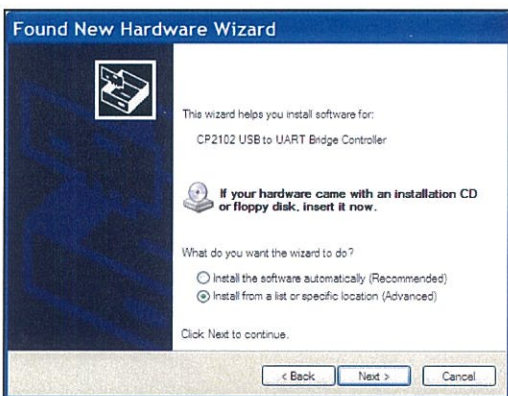
The window of "Found New Hardware Wizard" will show up again. Select "Install from a list or specific location". Click "Next".

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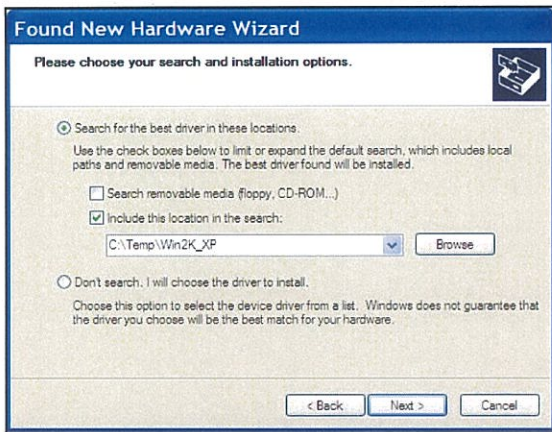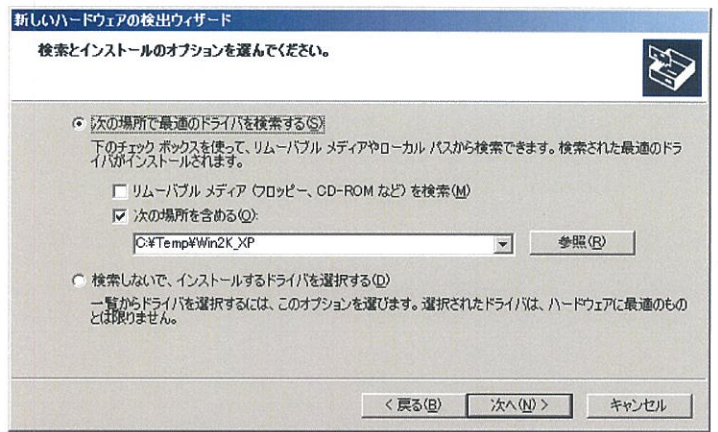
Select "Search for the best driver...".  
 Uncheck "Search removable media".  
 Check "Include this location in the search".  
 Click "Browse" and select "Win2K\_XP" folder.  
 Click "OK".  
 Click "Next".

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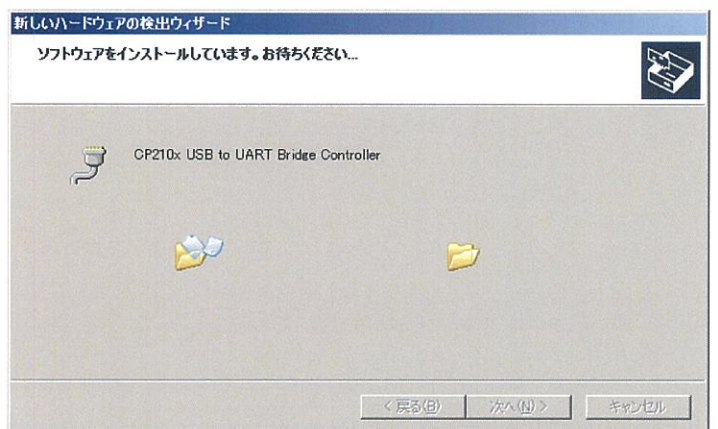
"CP210x USB to UART Bridge Controller" installation begins.

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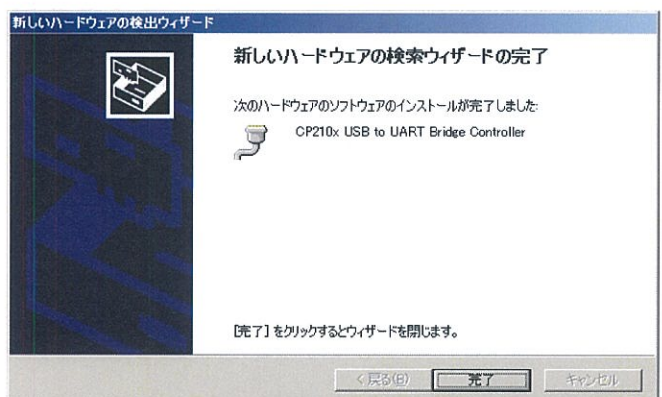
"CP210x USB to UART Bridge Controller" installation finished.  
 Click "Finish".

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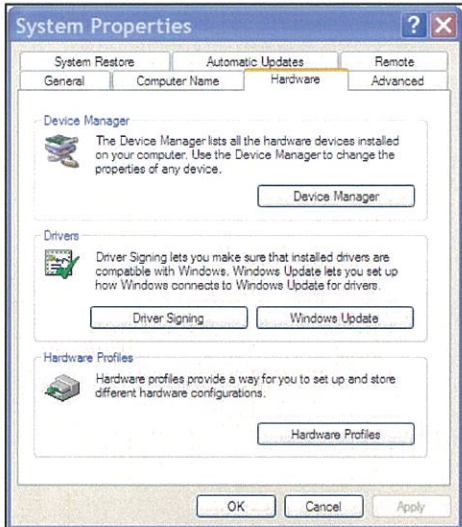
After connecting USB I/F cable and opening "Device Manager".


Go to "Start"→"Control Panel"→"System".  
Click "Hardware" tab.  
Click "Device Manager".

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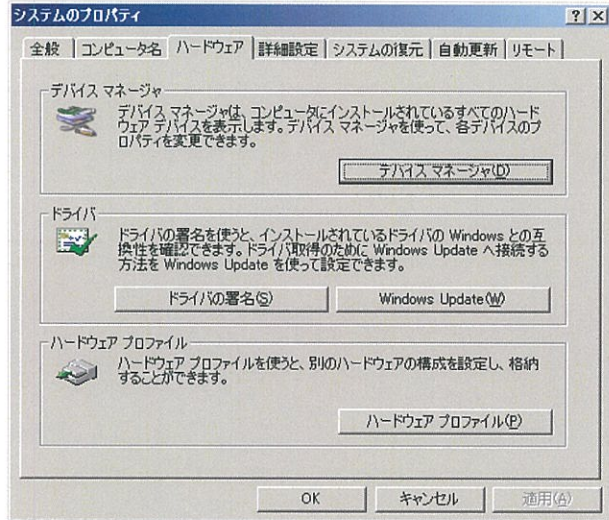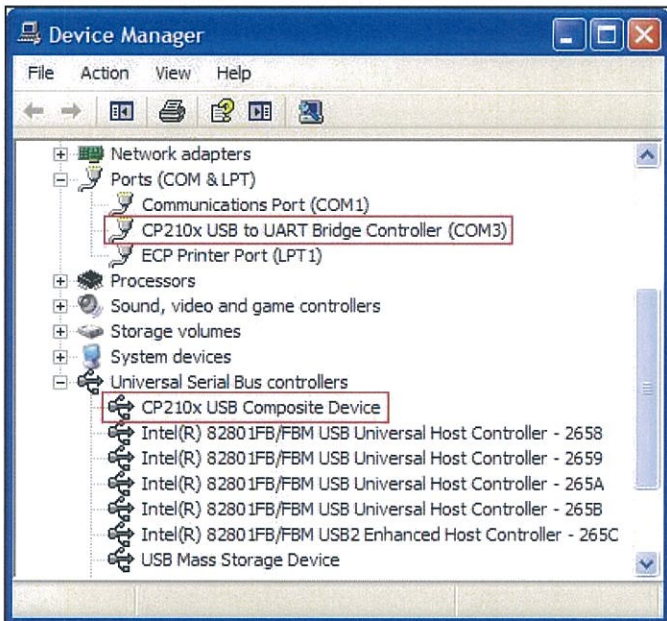
Confirm the port number and indication of "CP210x USB Composite Device" and "CP210x USB to UART Bridge Controller (COM\*)" as the window image.

This window shows that COM3 has the partition.

In EM Pro menu, go to "Set"→"Communication" then change the port to designated COM port.

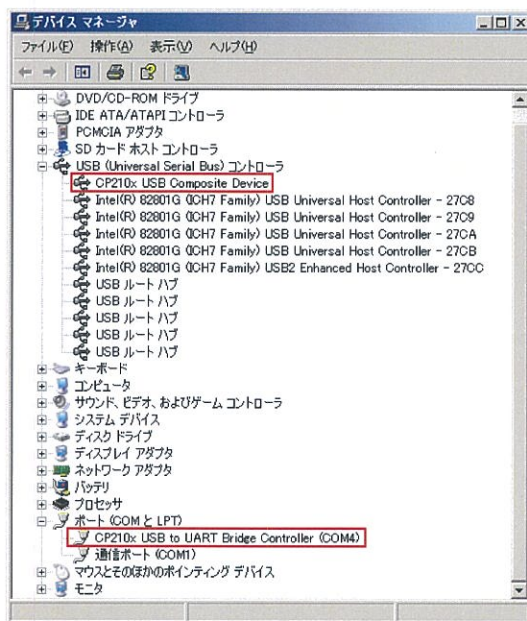After installation, the recommendation window of the driver update would show up, but ignore it.

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

## USB Driver uninstallation procedure

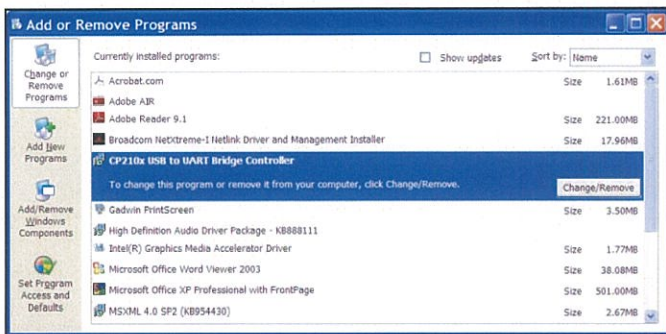
Go to "Start"→"Control Panel"→"Add or Remove programs".

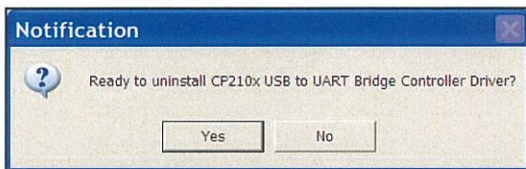Select "CP210x USB UART Bridge Controller".  
Click "Change/Remove".

Click "Yes".

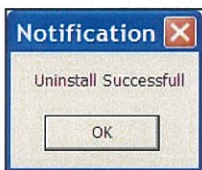Click "OK".

### CAUTION!

Specifications are subject to be changed without notice.

All trademarks used in this user manual are the property of their respective owners.

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Copyright Yoshimura Japan Co.,Ltd.
[Published by] YOSHIMURA JAPAN Co.,Ltd.